



BODY PLATE



CAPPING PLATE



FRAME PLATE



TAMPER



MIDDLE PLATE


POWDER SPREADER



SPILL GUARD



CAPSULE BODY & CAP

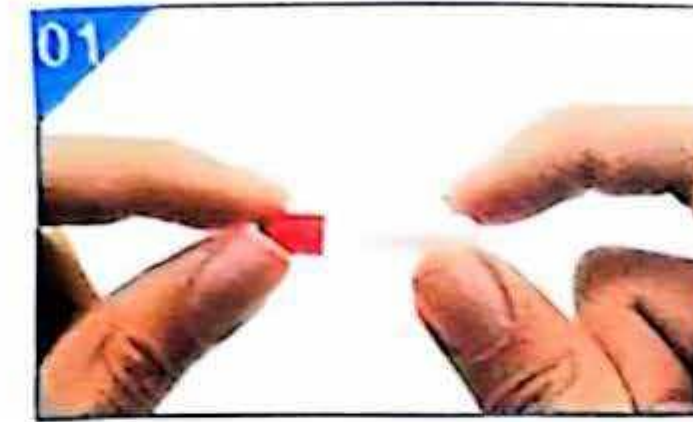

OPEN UP 100PCS OF YOUR CAPSULES AND PUT THEM IN SEPARATE BOWLS
The large part is the body and the small one is the cap



PLACE THE BODY PLATE ON A TABLE



PLACE THE FRAME PLATE ON TOP OF IT



POUR 100PCS OF THE CAPSULE BODY PARTS ON THE FRAME PLATE



SHAKE THE CPASULE MACHINE GENTLY
So that all the body parts of the capsule fall into their places.
If the capsule has fallen the wrong way around turn it by hand so all the capsule openings are facing up



LIFT UP THE FRAME PLATE AND PLACE THE SPILL GUARD




POUR YOUR POWDER ON TOP
Use the POWDER SPREADER to spread the powder evenly



USE THE TAMPER TOOL TO TAMP DOWN YOUR POWDER



ADD A BIT MORE OF YOUR POWDER AND FILL UP THE CAPSULES

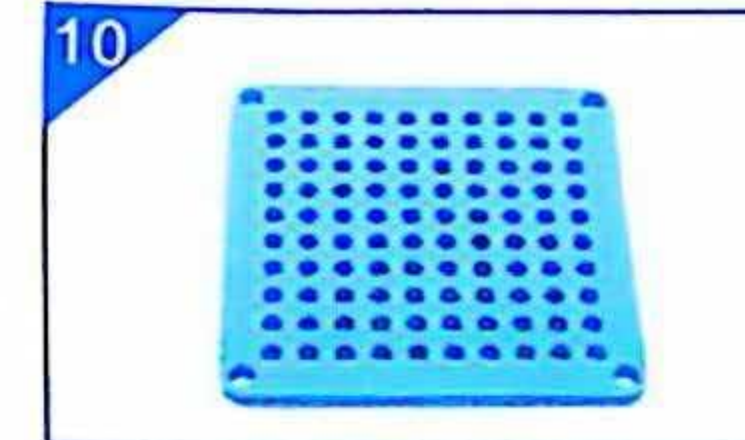


MOVE THE BODY PLATE A SIDE AND TAKE OUT THE CAPPING PLATE




PLACE THE FRAME PLATE ON TOP AND POUR 100PCS OF THE CAPSULE CAPS



SHAKE THE CAPPING PLATE GENTLY
So that all the capsule caps fall into their places. If the capsule has fallen the wrong way around turn it by hand so all the capsule openings are facing up

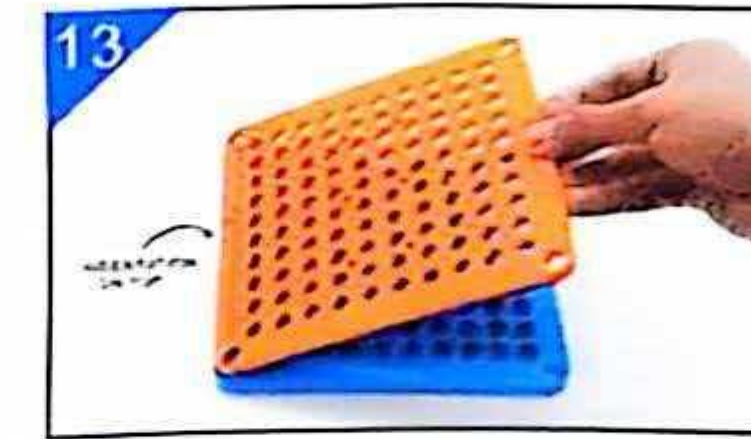


REMOVE THE FRAME PLATE AND PLACE THE MIDDLE PLATE ON TOP OF THE CAPPING PLATE

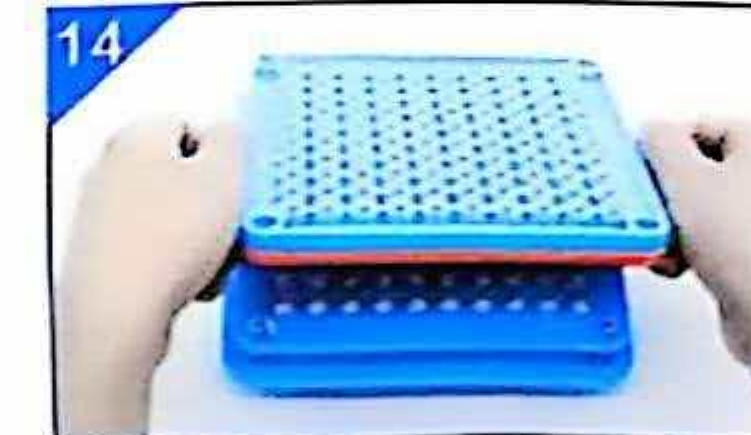


TAKE OUT THE BODY PLATE
FLIP THE CAPPING+MIDDLE PLATE ON TOP OF THE BODY PLATE



PRESS DOWN FIRMLY A COUPLE OF TIMES
Remember to press evenly all around the capsule machine.
You should hear multiple snapping sounds as the capsules close

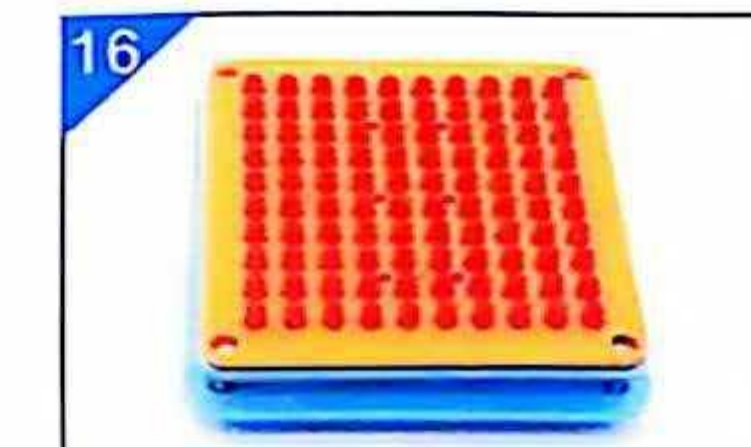

LIFT UP THE CAPPING PLATE

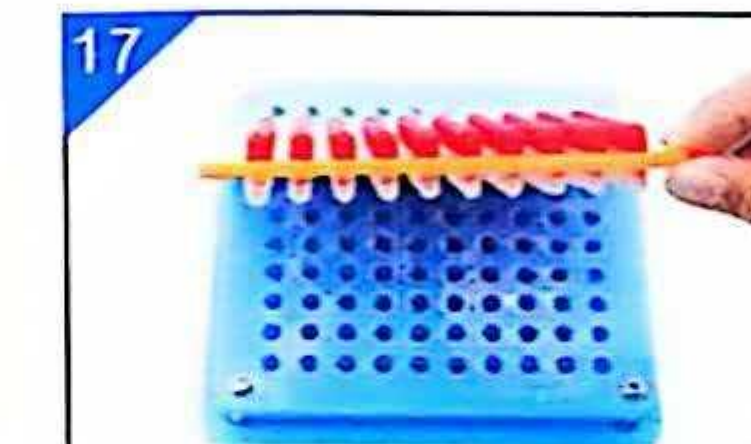

PICK UP THE MIDDLE PLATE AND FINISH YOUR CAPSULES!



PERFECT!